

Product Insert VARIANT Plex™-HT Lymphoma panel

VARIANT*Plex*-HT Lymphoma

Description

The VARIANTPlex-HT Lymphoma panel is a balanced pool of gene-specific primer (GSP) oligonucleotides that is optimized for use with VARIANT*Plex* reagents and molecular barcode (MBC) adapters to produce targeted NGS libraries. This product insert should be used in conjunction with VARIANT*Plex*-HT protocol for Illumina® (RA-DOC-058).

VARIANTPlex-HT Lymphoma contains **1787** GSPs targeting **51** genes commonly mutated in T- and B- cell lymphomas.

Description	Part number	Storage
VARIANTPlex-HT Lymphoma GSP1, 24 reactions or VARIANTPlex-HT Lymphoma GSP1, 96 reactions	SA24343241 or SA24343961	-20°C ± 10°C
VARIANTPlex-HT Lymphoma GSP2, 24 reactions or VARIANTPlex-HT Lymphoma GSP2, 96 reactions	SA24343242 or SA24343962	

Required reagent volumes

Protocol reference	Protocol step	Reagent	Volume per reaction (μL)
Α	Ligation Step 2 Elution	5mM NaOH	20
В	First PCR	VariantPlex-HT Lymphoma GSP1	8
С	First PCR	10mM Tris-HCl pH 8.0	18
D	First PCR	Purified PCR1 eluate	16
Е	Second PCR	VariantPlex-HT Lymphoma GSP2	8

2425 55th Street, Boulder, CO 80301 | archer-tech@idtdna.com

RA-DOC-067/REV02



VARIANT*Plex*™-HT Lymphoma panel

Recommended PCR cycling

	Step	Temperature (°C)	Time	Cycles
	1	95	3 min	1
	2	95	30 sec	
First PCR	3	60	10 sec	15
reaction	4	63	5 min (100% ramp rate)	_
	5	72	3 min	1
	6	4	Hold	1
	1	95	3 min	1
Second PCR reaction	2	95	30 sec	
	3	60	10 sec	20†
	4	65	5 min (100% ramp rate)	_
	5	72	3 min	1
	6	4	Hold	1

[†]The number of PCR2 cycles may be decreased if you regularly experience library yields greater than 200 nM.

Recommended reads and multiplexing

VARIANTPlex-HT Lymphoma libraries should be sequenced to a minimum of **5M reads**. Starting read depth recommendations for standard profiling may be adjusted based on user needs.

Archer™ Analysis settings

Sequencing data should be processed using Archer Analysis (v7.3 or greater). The VARIANTPlex-HT Lymphoma panel requires selection of the *SNV/Indel, Structural Variation*, and Copy Number Variation pipelines, found under the *DNA* Input Type (see the software user manual for further details on setting up analyses). Selection of the DNA Target Coverage pipeline is optional.

Processing of VARIANTPlex-HT Lymphoma libraries requires a one-time upload of the Custom Panel GTF. When performing DNA Target Coverage analysis, users must also select a Region

2425 55th Street, Boulder, CO 80301 | archer-tech@idtdna.com

RA-DOC-067/REV02



VARIANT*Plex*™-HT Lymphoma panel

of Interest BED file. Users may optionally add a Targeted Mutations VCF file for targeted SNV/Indel detection. Files can be obtained by contacting archer-tech@idtdna.com

Assay targets

Accession	Exon
NM_006015	1,2,3,4,5,6,7,8,9,10,11,12,13,14,15,16,17,18,19,20
NM_000051	1,2,3,5,6,7,8,9,10,11,12,13,14,15,16,17,18,19,20,21,22,23,24,25,26,27,28,29,30, 31,32,33,34,35,36,37,38,39,40,41,42,43,44,45,46,47,48,49,50,51,52,53,54,55,56, 57,58,59,60,61,62,63
NM_001351835	4
NM_004048	1,2,3
NM_000633	3
NM_000657	2
NM_001165	2,3,4,5,6,7,8,9
NM_004333	15
NM_000061	15
NM_032415	2,3,5,6,7,8,9,10,11,12,13,14,15,16,17,18,19,20,21,22,23,24,25
NM_053056	1,2,3,4,5
NM_006139	4
NM_006139	1,2,3,4
NM_001783	1,2,3,4,5
NM_000626	1,2,3,4,5,6
NM_000077	2,3
NM_001195132	3
NM_058195	1
NM_058197	1
NM_004936	2
	NM_006015 NM_000051 NM_001351835 NM_004048 NM_000633 NM_000657 NM_001165 NM_004333 NM_000061 NM_032415 NM_053056 NM_053056 NM_006139 NM_006139 NM_001783 NM_00077 NM_001195132 NM_058195 NM_058197

2425 55th Street, Boulder, CO 80301 | archer-tech@idtdna.com

RA-DOC-067/REV02



VARIANT*Plex*™-HT Lymphoma panel

Gene	Accession	Exon
CDKN2B	NM_078487	1
CREBBP	NM_004380	1,2,3,4,5,6,7,8,9,10,11,12,13,14,15,16,17,18,19,20,21,22,23,24,25,26,27,28,29, 30,31
CXCR4	NM_001008540	1
CXCR4	NM_001348056	2
CXCR4	NM_003467	1
DNMT3A	NM_001320893	1
DNMT3A	NM_022552	2,3,5,6,7,8,9,10,11,12,13,14,15,16,17,18,19,20,21,22,23
DNMT3A	NM_153759	1,2
DNMT3A	NM_175630	4
EP300	NM_001429	1,2,3,4,5,6,7,8,9,10,11,12,13,14,15,16,17,18,19,20,21,22,23,24,25,26,27,28,29, 30,31
EZH2	NM_004456	2,3,4,5,6,7,8,9,10,11,12,13,14,15,16,17,18,19,20
FOXO1	NM_002015	1,2
ID3	NM_002167	1,2
IDH2	NM_002168	4
JAK1	NM_002227	14,17,20,25
JAK3	NM_000215	3,10,11,12,13,14,15,16,17,18,19
KLF2	NM_016270	1,2,3
KMT2D	NM_003482	2,3,4,5,6,7,8,9,10,11,12,13,14,15,16,17,18,19,20,21,22,23,24,25,26,27,28,29,30, 31,32,33,34,35,36,37,38,39,40,41,42,43,44,45,46,47,48,49,50,51,52,53,54,55
MEF2B	NM_001145785	8
MEF2B	NM_005919	4,5,6,7,8,9,10
MYC	NM_002467	1,2,3
MYD88	NM_001172567	3
MYD88	NM_002468	4,5

2425 55th Street, Boulder, CO 80301 | archer-tech@idtdna.com

RA-DOC-067/REV02



VARIANT*Plex*™-HT Lymphoma panel

Gene	Accession	Exon
NFKBIE	NM_004556	1,2,3,4,5,6
NOTCH1	NM_017617	34, c.*370 to c.*380
NOTCH2	NM_024408	34
NRAS	NM_002524	2,3,4,5
PIM1	NM_001243186	1
PIM1	NM_002648	2,3,4,5,6
PLCG1	NM_002660	1,2,3,4,5,6,7,8,9,10,11,12,13,14,15,16,17,18,19,20,21,22,23,24,25,26,27,28,29, 30,31,32
PLCG2	NM_002661	2,3,4,5,6,7,8,9,10,11,12,13,14,15,16,17,18,19,20,21,22,23,24,25,26,27,28,29,30, 31,32,33
PTEN	NM_000314	1,2,3,4,5,6,7,8,9
PTEN	NM_001304717	1,2
PTPRD	NM_001171025	18
PTPRD	NM_002839	12,13,14,15,16,17,18,19,20,21,22,23,24,25,26,27,28,30,31,32,33,34,35,36,37,38, 39,40,41,42,43,44,45,46
PTPRD	NM_130392	18
PTPRD	NM_130393	12
RHOA	NM_001664	2,3
SF3B1	NM_012433	13,14,15,16,17,18,19,20,21
SOCS1	NM_003745	2
STAT3	NM_003150	19,20
STAT3	NM_139276	21
STAT5B	NM_012448	14,15,16,17
STAT6	NM_003153	8,9,10,11,12,13,14
TCF3	NM_001136139	18
TCF3	NM_003200	2,3,4,5,6,7,8,9,10,11,12,13,14,15,16,17,18,19

2425 55th Street, Boulder, CO 80301 | archer-tech@idtdna.com

RA-DOC-067/REV02



VARIANT*Plex*™-HT Lymphoma panel

Gene	Accession	Exon
TET2	NM_001127208	4,5,6,7,8,9,10,11
TET2	NM_017628	3
TNFAIP3	NM_006290	2,3,4,5,6,7,8,9
TNFRSF14	NM_003820	1,2,3,4,5,6,7,8
TP53	NM_000546	1,2,3,4,5,6,7,8,9,10
TP53	NM_001126113	10
TP53	NM_001126114	10
TRAF2	NM_021138	2,3,4,5,6,7,8,9,10,11
XPO1	NM_003400	15,16,18

Genes targeted for CNV

ARID1A	CD28	DNMT3A	KLF2	PIM1	SOCS1	TRAF
ATM	CD79A	EP300	KMT2D	PLCG1	STAT6	
B2M	CD79B	EZH2	MEF2B	PLCG2	TCF3	
BCL2	CDKN2A	FOXO1	MYC	PTEN	TET2	
BIRC3	CDKN2B	ID3	MYD88	PTPRD	TNFAIP3	
CARD11	CREBBP	JAK1	NFKBIE	RHOA	TNFRSF14	
CCND1	CXCR4	JAK3	NRAS	SF3B1	TP53	

Please contact archer-tech@idtdna.com to inquire about enabling additional genes for CNV detection.

SNPs and sites targeted for sample tracking

	rs560681	rs430046	rs987640	rs10776839	rs12393891
	rs740598	rs8078417	rs6444724	rs6530357	chrX 4429309
	rs1498553	rs9951171	rs6811238	rs5971553	chrX 11314433
	rs10773760	rs576261	rs13182883	rs5953060	chrY 6738552
	rs1058083	rs1109037	rs214955	rs6524626	chrY 19490214
	rs4530059	rs1523537	rs321198	rs5940270	
ľ	rs1821380	rs221956	rs4606077	rs722847	

SNPs may be used in combination to uniquely tag and track samples over time. Contact archer-tech@idtdna.com for further details.

2425 55th Street, Boulder, CO 80301 | archer-tech@idtdna.com

RA-DOC-067/REV02



Product Insert VARIANT*Plex*™-HT Lymphoma panel

Limitations of use

For research use only. Not for use in diagnostic procedures. Unless otherwise agreed to in writing, IDT does not intend these products to be used in clinical applications and does not warrant their fitness or suitability for any clinical diagnostic use. Purchaser is solely responsible for all decisions regarding the use of these products and any associated regulatory or legal obligations.

Safety data sheets pertaining to this product are available upon request.

This product or service is licensed under one or more of the following U.S. Patents: 8,835,358; 9,290,808; 9,290,809; 9,315,857; 9,708,659; and 9,816,137 owned by BD and is licensed solely for the use described in the associated product literature. No other rights, implied or otherwise, are granted to purchaser hereunder. Purchaser agrees, by way of example and not limitation, not to use this product to trace back the origin of a nucleic acid to an individual cell as a discrete entity (e.g., single cell analysis).

© 2023 Integrated DNA Technologies, Inc. All rights reserved. FUSIONPlex, VARIANTPlex, LIQUIDPlex, IMMUNOVerse, Archer Analysis, and Archer Assay Marketplace are trademarks of Integrated DNA Technologies, Inc. All other marks are the property of their respective owners. For specific trademark and licensing information, see www.idtdna.com/trademarks.

Revision History

Document Number	Date	Description of change
RA-DOC-067/REV01	July 2024	Initial release.
RA-DOC-067/REV02		Updated panel specification information to include PTEN and NRAS gene targets

2425 55th Street, Boulder, CO 80301 | archer-tech@idtdna.com

RA-DOC-067/REV02